



CATALYST TIMER RESET

ZAC-GC & ZAC-ULT Series of Zero Air Generators (Not ZAC-GC1500 Zero Air Generators)

The catalyst has a limited lifespan and needs replacement after three years of operation (please see the operations manual). As a reminder to the operator, the generator will display a warning message “Catalyst life alarm”, when the life of the catalyst is exceeded. The red external “check system” LED begins to flash. From this time on, the operator has 60 days to replace the catalyst or the control board disables the system.

After replacement of the catalyst, the timer needs to be reset.

- 1) Turn off the power to the unit.
- 2) Locate the jumper (JP1) just on the right side of the display (DS1) on the bottom left hand side of the main circuit board.
- 3) Move the jumper from the (Pin-1 / Pin-2) position to the (Pin-2 / Pin-3) position.
- 4) Turn on the power to the unit and leave on for 30 seconds
- 5) Turn off the power to the unit.
- 6) Move the jumper from the (Pin-2 / Pin-3) position to the (Pin-1 / Pin-2) position.
- 7) Turn on the power to the unit. The catalyst timer has been reset.

